

NOMAD NANO

High Flow Pulsatile Pump System

A powerful, compact, precision-engineered pulsatile flow system built for R&D, engineering, and field demonstrations. The NOMAD NANO is the smallest and most economical pulsatile pump in the market, delivering flow rates up to 8 L/min while maintaining the performance, versatility, and control required for cardio and neurovascular simulation.




System Features

- Pulse Rate: 0-180 BPM (250ms pulse duration) • Adjustable Flow Power (0-100% or 0-5L/min) • Can Run Continuous
- Maximum Output Flow: 8L/min (1/2" Barbed Ports) • Maximum Output Pressure: ~450 mmHg • Travel-Sized System
- Configurable Ports: Valved (CPC) or Barbed Ports • Low Power Rating 12V System (Power Supply: 12VDC 5A - 60W)

System Specifications

- **Dimensions:** (5.00*4.00*5.00)in • **Enclosure Materials:** ABS or PLA • **System Total Weight:** 1.5lb


Flow and Pressure Waveform Characterization Test

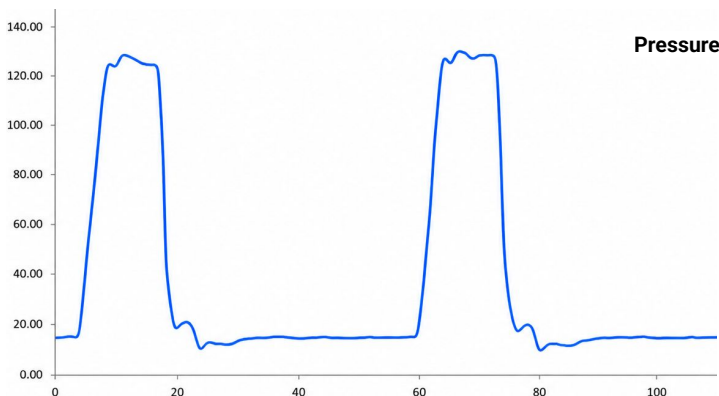

Test Configuration

Pressure and flow testing was performed with the NOMAD NANO connected in a closed-loop configuration using a 12-inch, 3/8-inch ID rigid PVC circuit. A water reservoir was filled with one liter of water, thermally conditioned to 98.6°F (37°C). The system was operated at 60 BPM and 100% Flow Rate (CPC - PLCD 16006).

Pressure waveform data was acquired using the Vascular-Labs DPM (Digital Pressure Monitor). Flow rate data was collected using a multi-sensor array comprising an analog flow meter, a digital flow sensor, and a high-speed camera.

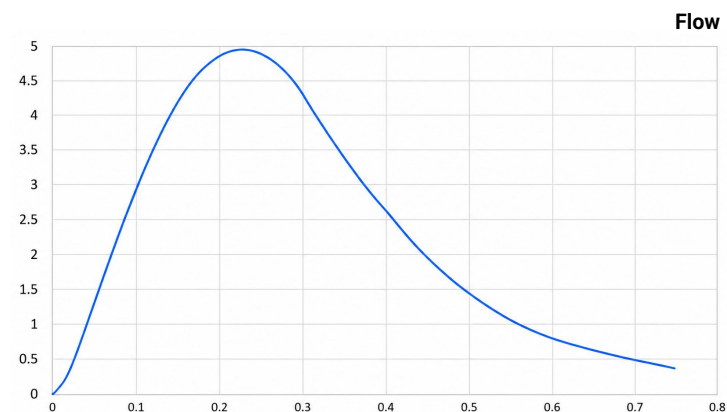

Disclaimer: The content of this document is believed to be correct at the time of release. However, Vascular-Labs and its affiliates offer no guarantees whatsoever, express or implied, in case of typographic or other errors. All specifications are subject to change without notice. Vascular-Labs products are designed for training purposes and are not classified as medical devices. They are not intended for diagnosing treating, curing, or preventing any medical condition, and their safety and efficacy as medical devices have not been established.